




Classic Stills Accessories



Floor Stand for Cylindrical Storage Tank

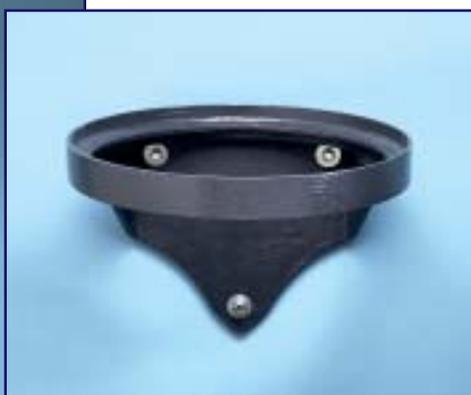

Wall Bracket for Cylindrical Storage Tank

Floor or Wall Mounted Storage!

- Mount storage tanks on a wall or place them on a stand
- Maintain sterility of stored water with a UV lamp

Ultraviolet Lamp Assembly

- The sterility in the tank is maintained with the UV lamp.
- The ultraviolet assembly consists of a cylindrical electrical housing which includes a starter, ON/OFF switch, 6 foot electrical cord, lamp and ballast.
- The housing is designed to be mounted on top of a storage tank with the lamp extended vertically into the tank.
- The lamp is enclosed by a tubular, transparent, inert sheath. The sheath is sealed at the neck by a grommet installed on top of the storage tank. When it is necessary to change the lamp, only the lamp need be removed. The reservoir cover and sheath remain in place at all times so distilled water is never exposed to possible contamination from the air.
- The ultraviolet lamp operates from 120 VAC, 50/60 Hz, 35 Watt electrical power.

Each ultraviolet lamp assembly is furnished with one bulb which should be replaced every 90 days if operated continuously. Rectangular tanks are always provided with Ultraviolet Assembly.

Product Description

Floor Stands for 10-50 Gallon Cylindrical Storage Tanks

- Floor stands for 10-50 gallon storage tanks are fabricated from heavy gauge sheet metal and are provided with adjustable feet.
- Floor stands for 100-200 gallon storage tanks are fabricated of thick steel angle-iron.

Wall Bracket for Cylindrical Storage Tanks

- Wall brackets for 10 gallon tanks are fabricated from heavy gauge aluminum.
- Wall brackets for 25 gallon tanks are fabricated of thick steel angle-iron.

ORDERING INFORMATION

Floor Stands for Cylindrical Storage Tanks					
Model #	Size of tanks to be held (gallon)	Overall Dimensions			US List Price
		W	H	D	
H1000	10	11 (28)	27 (69)	11 (28)	\$458.00
H1001	25	19 (48)	27 (69)	19 (48)	\$493.00
H1002	50	22 (56)	20 (51)	22 (56)	\$557.00
H1003	100	29 (74)	11 (28)	29 (74)	\$680.00
H3230	200	39 (99)	22 (56)	39 (99)	\$1,079.00
Ultraviolet and Lamp Assembly					
Model #	Description				US List Price
H3240	Wall bracket for cylindrical storage tanks (10 gal)				\$414.00
H3242	Wall bracket for cylindrical storage tanks (25 gal)				\$673.00
H4005	UV Lamp Assembly and Lamp				\$729.00
04141	Replacement Ultraviolet Lamp for 120V operation				\$54.00